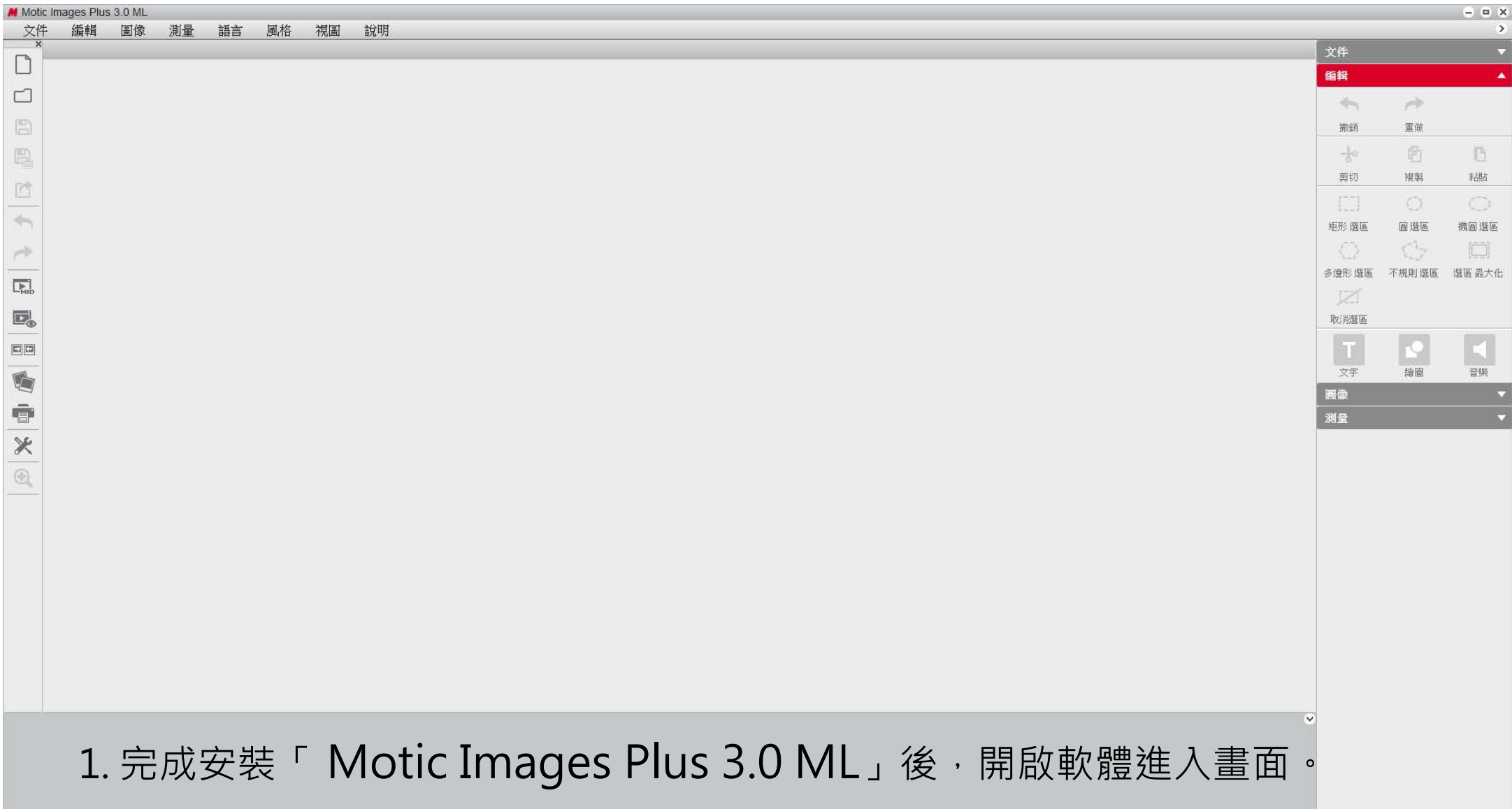
Motic[®]

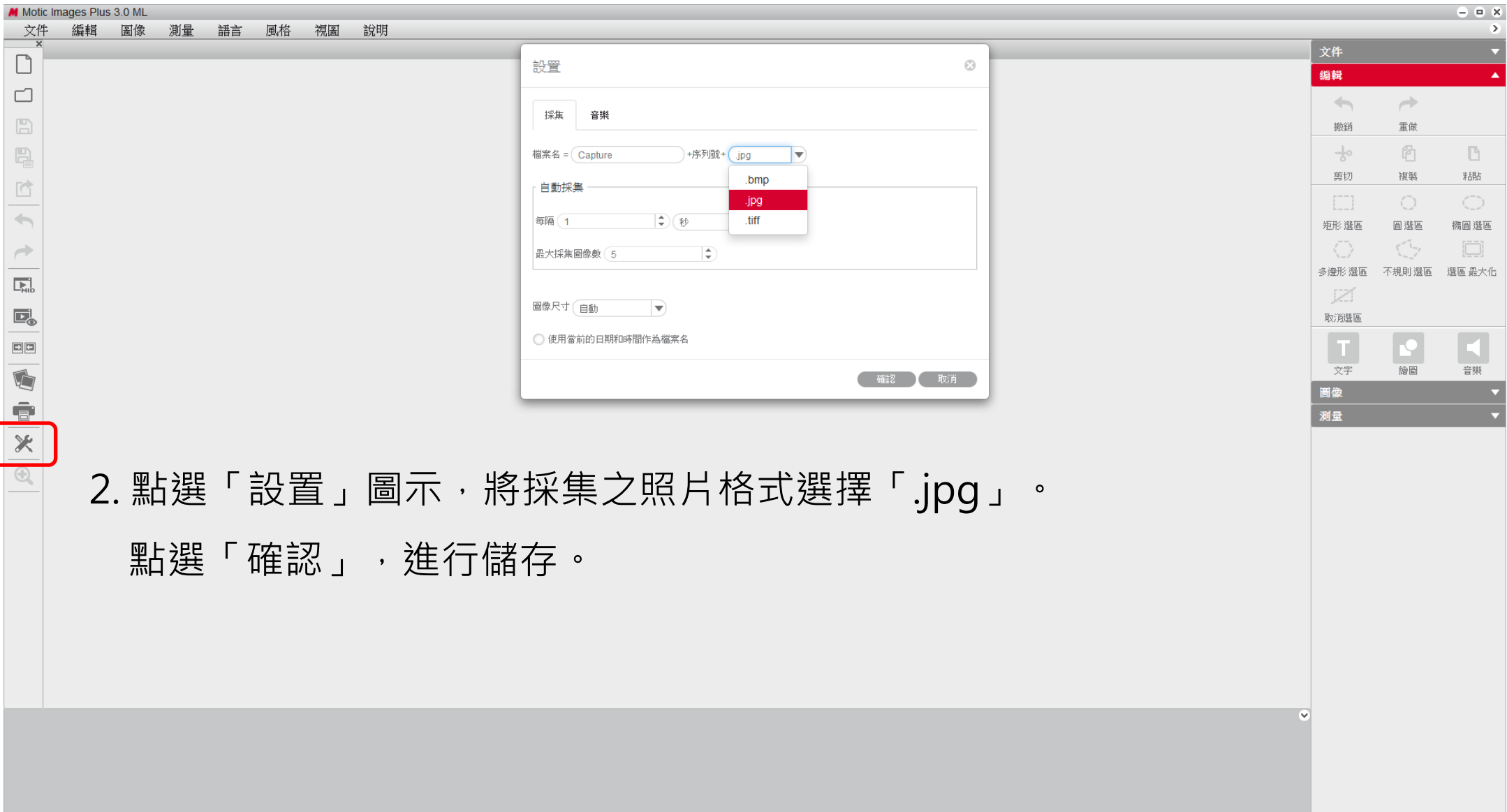
Motic Images Plus 3.0 ML

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The screenshot displays the Motic software interface. The main window title is 'Motic 動態成像模組'. The top menu bar includes '說明', '語言', and '風格'. Below the menu bar, there are controls for '信數' (100X), '顯示信數', and 'ROI 預覽'. The central area is a large gray rectangle. On the right side, there is a '基本調節' (Basic Adjustment) panel. The '視頻設備' (Video Device) dropdown menu is open, showing three options: 'Motic Demo', 'Motic 5.0MP' (highlighted in red), and 'Motic Wifi'. Other settings in the panel include '分辨率', '曝光', '增益', '偏移', '增強', '伽馬', and buttons for '設置', '計算白平衡', '讀取背景', '鏡像', '倒置', '充滿視窗', and '全屏'.

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注意：Motic Demo / Motic Wifi 為軟體預設裝置

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The screenshot displays the Motic software interface. The main window shows a dark circular field of view, likely representing a microscope's field of view. The interface includes a top menu bar with options like '說明', '語言', and '風格'. Below the menu bar, there are status indicators for 'FPS:8.91', '倍數: 100X', '顯示倍數', 'ROI 預覽', and 'Plus_100X(1, 1)'. On the right side, there is a '基本調節' (Basic Adjustment) panel with various settings:

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
The screenshot shows the Motic software interface. The main window displays a large blue circle on a dark background. A calibration window titled '校準' is open, showing a smaller blue circle and various settings. The '高級設置' (Advanced Settings) panel on the right is also visible.

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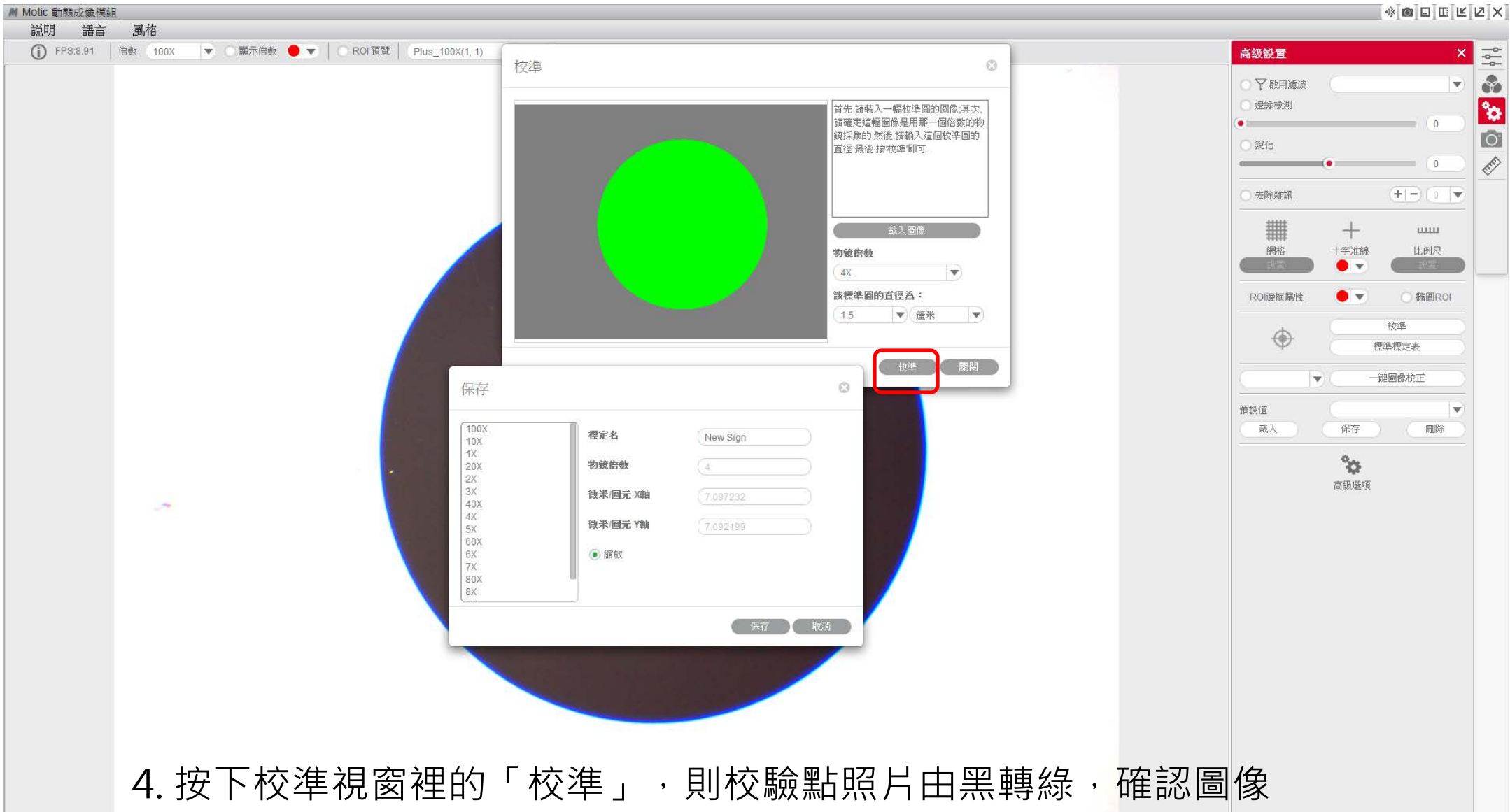The screenshot shows the Motic software interface with the 'Calibration' dialog box open. The dialog box contains a central image of a green circle on a black background. To the right of the image, there is a text box with instructions: '首先,請裝入一幅校準圖的圖像,其次,請確定這幅圖像是用那一個倍數的物鏡採集的,然後,請輸入這個校準圖的直徑,最後,按校準即可。' Below this text are fields for '載入圖像', '物鏡倍數' (set to 4X), and '該標準圖的直徑為:' (set to 1.5 厘米). At the bottom of the dialog, the '校準' button is highlighted with a red box. In the foreground, the '保存' dialog box is open, showing a list of magnification options (100X to 8X) and fields for '標定名' (New Sign), '物鏡倍數' (4), '微米/圖元 X軸' (7.097232), and '微米/圖元 Y軸' (7.092199). The '縮放' option is selected. The '保存' button is highlighted.

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